



KFU

جامعة الملك فيصل

KING FAISAL UNIVERSITY

جامعة ووطن.. نماء.. واستدامة..



INSTRUMENT STANDARD OPERATING PROCEDURE MANUAL

College of Medicine



SAFETY AND LABORATORY COOMMITTEE, C.O.M, K.F.U.

Prepared by

Document Number	Name	Signature	Date
CM/EMU-011	husain musa al nwaiser ali althani		26/4/2021
Revision Number	Approved by		
R0	Dean:		
Write the revision number	Department:		

Document History

Document Section	Details of Amendments	Date	Modified by (Initials)
SOP	First Draft on SOP for the operation of Leica EM SCD500 – High Vacuum Sputter Coater		

1. OBJECTIVE

- ☐ The document describes the operation of **Leica EM SCD500 – High Vacuum Sputter Coater for SEM**
 - Specimens for SEM is coated with sputtered metal atoms using argon gas as ionized and accelerated onto the target's surface.

2. SCOPE

- ☐ Full automated software processor system
- ☐ Process gas (argon is fed at 0.5 mbar pressure).
- ☐ Sputter working distance 50mm.
- ☐ After specimen processing must be coated using Platinum ,Silver or Gold or Carbon (layer thickness 10 nm).

3. RESPONSIBILITIES

- ☐ It is the responsibility of designated personnel in the lab to train staff and students on this procedure and to ensure adherence to this procedure under supervision.
- ☐ It is the responsibility of designated personnel (staff or Student) to follow the instructions of this procedure under supervision.

4. REFERENCES

- ☐ It is the responsibility of designated personnel in Research lab to train staff and students on this procedure and to ensure adherence to this procedure under supervision.
- ☐ It is the responsibility of designated personnel (staff or Student) to follow the instructions of this procedure under supervision.

5. DEFINITIONS

- ☐ Closed fully automated **Leica EM SCD500 – High Vacuum Sputter Coater**
- ☐ Controlled programming software.
- ☐ Housing with control, display and process selection panel
- ☐ Vacuum system(turbomolecular pump,roughing pump and vacuum ion gauge).
- ☐ Specimen chamber with build in height adjustable table.
- ☐ Gas inlet system with manually controlled gas dosing valve,electromagnetically controlled cut-off valve and push button activating rinsing gas.

- ☐ Hinged sputter head with planar magnetron magnet system, target holder and providing shield .
- ☐ HT 010 high voltage supply for sputter and etching device.
- ☐ Electromagnetically controlled venting valve coupled with main switch.
- ☐ Carbon tips(6.15,30mm long
- ☐ Platinum wire(0,1 mm or 0,2mm)
- ☐ Chrome coated tungsten wire.

6. SAFETY PRECAUTIONS

- ☐ Tightening the gas dosing valve with force will damage the valve seat.
- ☐ Before switching on the electricity check all the flanges are closed.
- ☐ It is necessary to switch on the **High Vacuum Sputter Coater EM SCD 500** after service work or after longer non- use.
- ☐ It is necessary to switch off the **High Vacuum Sputter Coater EM SCD 500** only for service operations before periods of long term non- use
- ☐ Water cooling is not required if sputter time is short (less than 300 seconds and sputter below 60Ma.
- ☐ Rinsing is not required if gas lines are filled with argon.
- ☐ Do not over tighten the gas dosing valve to avoid the damage.
- ☐ Do not adjust gas dosing valve while turbo pump is slowing down.

7. PROCEDURE FOR OPERATING Leica EM SCD500 – High Vacuum Sputter Coater

7.1. Turning on the instrument:

- ☐ Adjust gas cylinder
- ☐ Switch ON the System.

7.2. Software Operation.

7.3. Steps of the Procedure

- ☐ Evacuation of vacuum chamber.
- ☐ It is necessary to switch off the **High Vacuum Sputter Coater EM SCD 500** only for service operations before periods of long term non- use.
- ☐ Sample loading after adjust the sample table
- ☐ Install glass chamber and safety cover
- ☐ Close the vacuum chamber
- ☐ Switch on the system .
- ☐ Adjust the time and sputter current (100mA).
- ☐ Wait for vacuum.
- ☐ Sputtering with full thickness control
- ☐ Wait for vacuum.
- ☐ Switch off system after sputtering process.

- ❑ **Evacuate specimen chamber.**

7.4. Turning off the Instrument

END OF PROGRAM

- **To turn off the instrument.**
- Evacuate specimen chamber
- Click shut down of computer
- Turn off the software.
- Turn off the electric main power source.
- All panel displays go out.
- Turn off the main switch

7.5. Warning

Write any the warning in BOLD and RED FONT.

